

1 Weight sets

- Make sure that the case is tightly closed and / or fix the closure
- Ideally pack the case in a bag. This prevents weights from being lost if the closure should open during transport
- Place the case with the bag in a sufficiently large outer package
- Pad these with additional filling material to give the case a stable position

2 Single weights

- If necessary, pad out the weight in the case with additional filling material so that it is well protected during shocks
- Make sure that the case is tightly closed and / or fix the closure
- Ideally pack the case in a bag. This prevents weights from being lost if the closure should open during transport
- Place the case with the bag in a sufficiently large outer package
- Pad these with additional filling material to give the case a stable position

3 Balances and other measuring devices

- **In principle: always follow the manufacturer's instructions**
- Use the original packaging for shipping
- Remove the weighing plate and place it separately in the package
- If feasible, install the transport safety
- For analytical balances, fix the doors of the protective cover, so that they remain tightly closed during transport
- **Please send all parts required for operation (e.g. the power supply unit)**

Kindly enclose the order form and any additional documents (delivery note, previous calibration certificate etc.)

Kindly address the shipment to:

KERN & SOHN GmbH
Calibration laboratory
Ziegelei 1
72336 Balingen

The listed notes are only to be considered as recommendations and can not be transferred to all devices. Shipping is always at your own risk.